
MODE D'EMPLOI // INSTRUCTIONS FOR USE
DMC 2000S

FRANÇAIS :

ALLUMAGE DU DOSIMETRE :

Appuyer **une fois** sur le bouton de commande et relâcher,
→ Il apparaît successivement à l'écran : « **CHANGE** » puis « **ENTER** »
Lorsque ce message « **ENTER** » est à l'écran, **appuyer** immédiatement sur le bouton de commande.
→ Le dosimètre est allumé

MISE EN VEILLE DU DOSIMETRE :

Appuyer **longuement**, sans relâcher, sur le bouton de commande,
→ Il apparaît successivement à l'écran : « **CHANGE** » puis « **GO OUT** »
Lorsque ce message « **GO OUT** » est indiqué, **relâcher** immédiatement le bouton de commande.
→ Le message «**OUT**» apparaît brièvement et le DMC se met ensuite en pause.
Le nom d'utilisateur (ex. SPS) est affiché quand le dosimètre est en pause.

ENGLISH :

SWITCHING TO THE MEASUREMENT MODE

Press once the selector button and release it,
→ The following messages are displayed in succession: « **CHANGE** » then « **ENTER** »
While « **ENTER** » is displayed, **press** immediately the selector button,
→ The DMC is activated

SWITCHING TO THE PAUSE MODE

The selector button should be **held pressed** (don't release it),
→ The following messages are displayed in succession: « **CHANGE** » then
« **GO OUT** »
When this « **GO OUT** » message is displayed, immediately **release** the selector button,
→ The message « **OUT** » appears briefly on the display and the DMC is in the pause mode.
An assigned name (i.e. SPS) is displayed.